F-SAFE

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SST-... 12/20

SST-... 6/20

SST-... 11/6

DE – Montage- und Bedienungsanleitung	2
EN – Installation and Operating Manual	36
FR – Instructions de montage et d'emploi	70
GS-Zertifikate	103

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Installation and Operating Manual

Hazardous-substance Cabinets

Table of contents

1	Important notes	40
1.1	General safety-related notes	41
1.2	Intended use	41
1.3	Functional testing, safety-related review, care, and maintenance	41
1.4	Warranty, declaration of transfer	42
1.5	Property rights	42
2	Technical data	43
3	Transport	43
4	Set-up/Initial start-up	44
4.1	General	44
4.2	Aligning the products	44
4.3	Tilt safety device	45
4.4	Base panel	45
4.5	Removing the transport restraint	46
5	Closure types/Locks	47
6	Interior fittings	47
6.1	Base tray	47
6.2	Tray bases	48
6.3	Pull-outs	48
7	Storage	48
8	Earthing	49
9	Ventilation	50
10	In case of fire	50
10.1	Opening the product after the fire	50
11	Disposal	50
12	Connection diagrams	51
13	Optional accessories	52
13.1	General safety information	52
13.2	Intended use	52
13.3	Warranty, surrender of goods	53
13.4	Single-cable / single-pipe penetration	54
13.5	Fire-resistant cable / pipe bulkhead	56
13.6	Recirculation fan, type SST-UL V4	58
13.6.1	Instructions for safe use	58
13.6.2	Specific hazards	59
13.6.3	Operating principle, safe installation, commissioning, adjustment, installation / removal, maintenance	59

13.6.4	Troubleshooting	61
13.6.5	Specifications, characteristics	61
13.6.6	Accessories shipped with the SST-UL	63
13.6.7	Installing the SST-UL.....	63
13.7	Exhaust fan, type SST AL	64
13.7.1	Instructions for safe use	64
13.7.2	Specific hazards	65
13.7.3	Safe installation, commissioning, adjustment, installation / removal, maintenance	65
13.7.4	Troubleshooting	66
13.7.5	Specifications.....	66
13.7.6	Accessories shipped with the SST-AL	68
13.7.7	Installing the SST-AL for cabinets	68

Dear customer,

You've decided in favour of a high-quality PROTECTO product.

Thank you very much.

Please carefully observe the operating manual below so we can guarantee that your whole system functions perfectly.

We can't undertake any warranty with non-observance.

We expressly reserve the right—even without separate notice—to make technical changes that serve to improve our products or are occasioned by legal changes.

This operation/installation manual may not be reprinted or reproduced—even in part—without the written consent of PROTECTOPLUS GmbH.

PROTECTOPLUS GmbH, Albert-Betz-Straße 1, 24783 Osterrönfeld

Tel.: +49 4331 4516-0 | Fax: +49 4331 4516-11

info@protecto.de | www.protecto.de

1 Important notes

Setting up and installing the product to be plumb and horizontal is imperative!





Individual parts must not be set down on their corners; storage area must be covered in advance with soft material such as a packing blanket!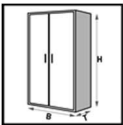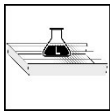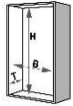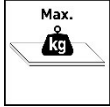

Caution! Hand injuries while actuating the pull-out may occur in the area of the guide rail.

Legend signs

	Exterior dimensions		Max. base-tray container size
	Interior dimensions		Max. lower-shelf load
	Total empty weight		Volume and pressure loss during ten air changes
	Max. mass		Safety cabinet under DIN EN 14470-1
	Max. base-tray collection volume		

1.1 General safety-related notes

- The operating/installation manual must be constantly stored where the product is being used.
- Legally prescribed safety notices are to be attached to places visible to the users.
- All safety and hazard notices, as well as the name plate, must be kept in legible condition.
- Use the product only when it's in proper condition.
- Observe the technical data in our catalogue or data sheet for each product.
- Improper use can impair the protection function.
- The products must not be stored outdoors and must be protected from water, penetrating moisture or splashing water. The products are not suitable for damp rooms.
- Observe all of the instructions in this operating and installation manual.
- Listing all of the valid regulations and directives would exceed the scope of this document. So always check to see whether your system meets valid regulations.

1.2 Intended use

- The products described in this operation/installation manual are built according to the latest state of technology and recognized safety-related rules.
- They may only be used:
 - as intended and
 - in flawless safety-related condition.
- Any other use is considered improper. PROTECTO assumes no liability for damage and the resulting consequential damage arising from this.

1.3 Functional testing, safety-related review, care, and maintenance

PROTECTO safety cabinets are safety-related systems.

Pursuant to sec. 4, para. 3 of the Workplaces Ordinance, ArbStättV, sec. 10 of the Ordinance on Industrial Safety, and Trade Associations rule BGR 234, a technical expert must conduct at least one annual safety-related inspection.

PROTECTO offers both individual inspections as well as maintenance agreements.

Daily inspection:

- Liquids must be immediately collected and properly disposed of.

Monthly inspection:

- Locking systems (cylinder lock)
- Hinges
- Door closers (if present)
- Correct seating and condition of the fire-protection seals
- Inside the exhaust and ventilation openings

1.4 Warranty, declaration of transfer

The installation instructions must be strictly adhered to in order to guarantee our products' optimal functioning.

PROTECTO assumes no liability for damage to other facilities that may arise from the installation or use of the product.

The warranty from PROTECTO extends to the delivered products.

Changes or modifications to the design may be conducted only after consultation with PROTECTO; otherwise the certification/warranty will be invalidated.

The installer assumes the warranty for installation services.




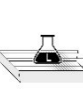
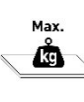
Warranty and liability claim for personal injury and property damage are excluded if they can be traced back to one or more of the following causes:

- Improper assembly, start-up, operation, and maintenance
- Instructions concerning transport, storage, operation, and assembly being disregarded
- Improper repairs
- The occurrence of emergencies caused by third parties' external actions, and force majeure.

1.5 Property rights

Utility models have been filed with the German Patent Office in order to safeguard innovation and design.

2 Technical data

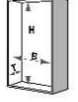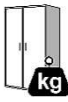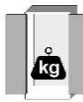
 Item	 HxWxD			 HxWxD			 About	 About	 About	 About	 About	 m ³ /h, Pa
	H	B	D	H	B	D	kg	kg	kg	kg	kg	10
6/20 F30	1960	595	595	1600	475	515	147	**	19	17	75	5 / About 1
6/20 F30 (4 p.-o.) (5 pull-outs) (6 pull-outs)	1960	595	595	1600	475	515	206	506	*	*	75	5 / About 1
							220	595				
							235	685				
12/20 F30	1960	1195	595	1600	1075	515	226	**	44	40	75	10 / About 1
12/20 F30 (4 p.-o.) (5 pull-outs) (6 pull-outs)	1960	1195	595	1600	1075	515	291	591	*	*	75	10 / About 1
							308	683				
							324	774				
11/6 F30	635	1100	550	375	980	472	97	**	24	21	75	5 / About 1
11/6 F30	635	1100	550	375	980	472	97	172	*	*	75	5 / About 1
6/20 F90	1960	600	595	1603	516	483	340	**	20	18	75	5 / About 1
6/20 F90 (4 p.-o.) (5 pull-outs) (6 pull-outs)	1960	600	595	1603	516	483	394	694	*	*	75	5 / About 1
							408	783				
							422	872				
12/20 F90	1960	1200	595	1603	1116	483	600	**	44	40	75	10 / About 1
12/20 F90 (4 p.-o.) (5 pull-outs) (6 pull-outs)	1960	1200	595	1603	1116	483	673	973	*	*	75	10 / About 1
							691	1066				
							710	1160				
11/6 F90	661	1100	500	491	1016	355	185	**	24	21	75	5 / About 1
1/6 F90	661	1100	500	491	1016	355	185	260	*	*	75	5 / About 1

* See the enclosed manual excerpts

** + max. 75 kg per storage level

Technical changes are reserved!


3 Transport

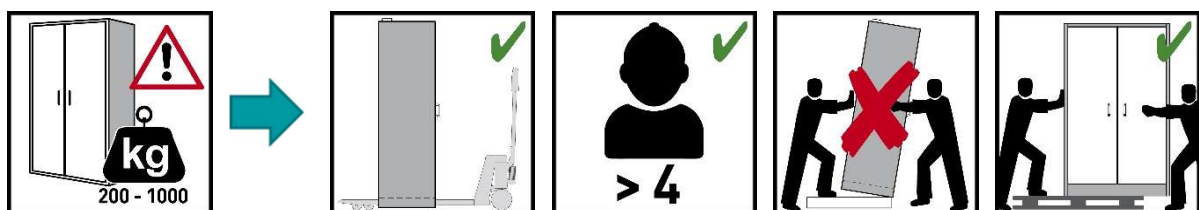

Use a forklift to transport the enclosure upright, tied up, and secured against slippage to the final set-up location. The transport restraints in the door joints may only be removed directly at the set-up location! The products are only functional after all of the transport restraints are removed. Improper transport can lead to hidden damage to the fire-protection insulation.

The cabinet may only be tilted smoothly and over its full area!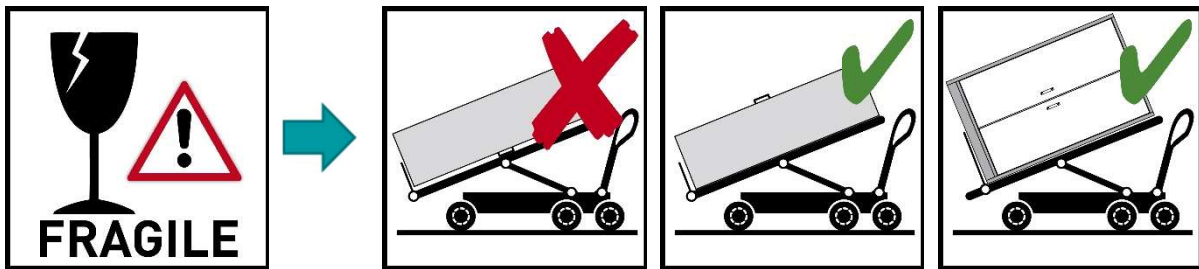



4 Set-up/Initial start-up

4.1 General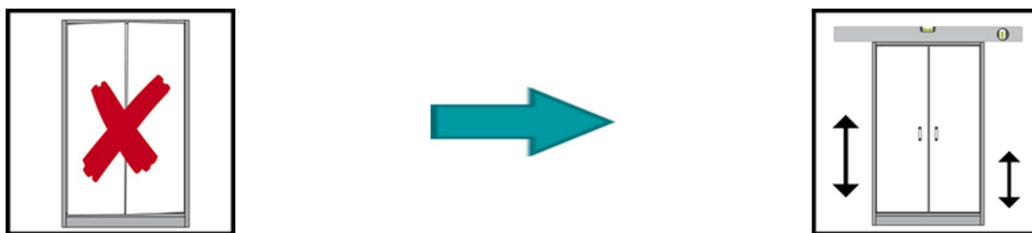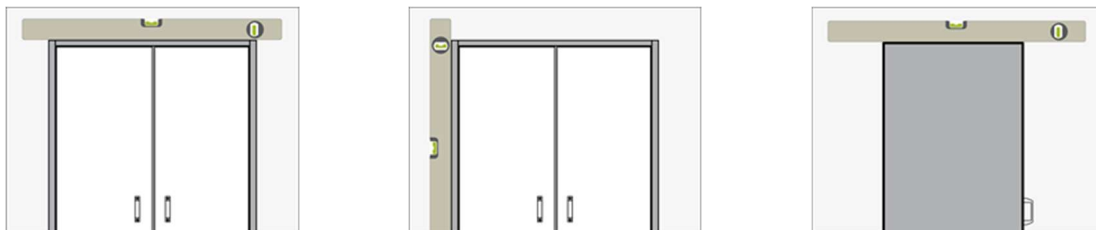

Technical changes are reserved!

4.2 Aligning the products
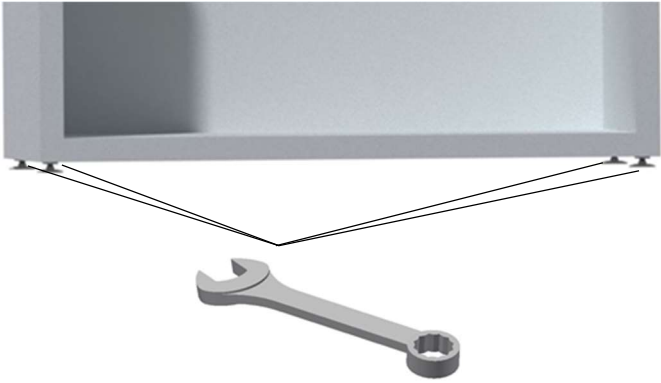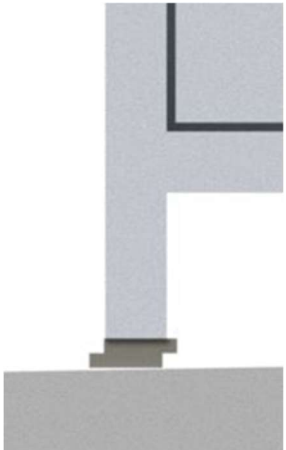
<p>6/20 F30 12/20 F30 6/20 F90 12/20 F90</p>		
<p>11/6 F90</p>		
<p>11/6 F30</p>		

4.3 Tilt safety device

Is needed only in products with pull-outs—please obtain information from the enclosed operating manual here.


Technical changes are reserved!

4.4 Base panel

Base and screws are included in the cabinet.

4.5 Removing the transport restraint

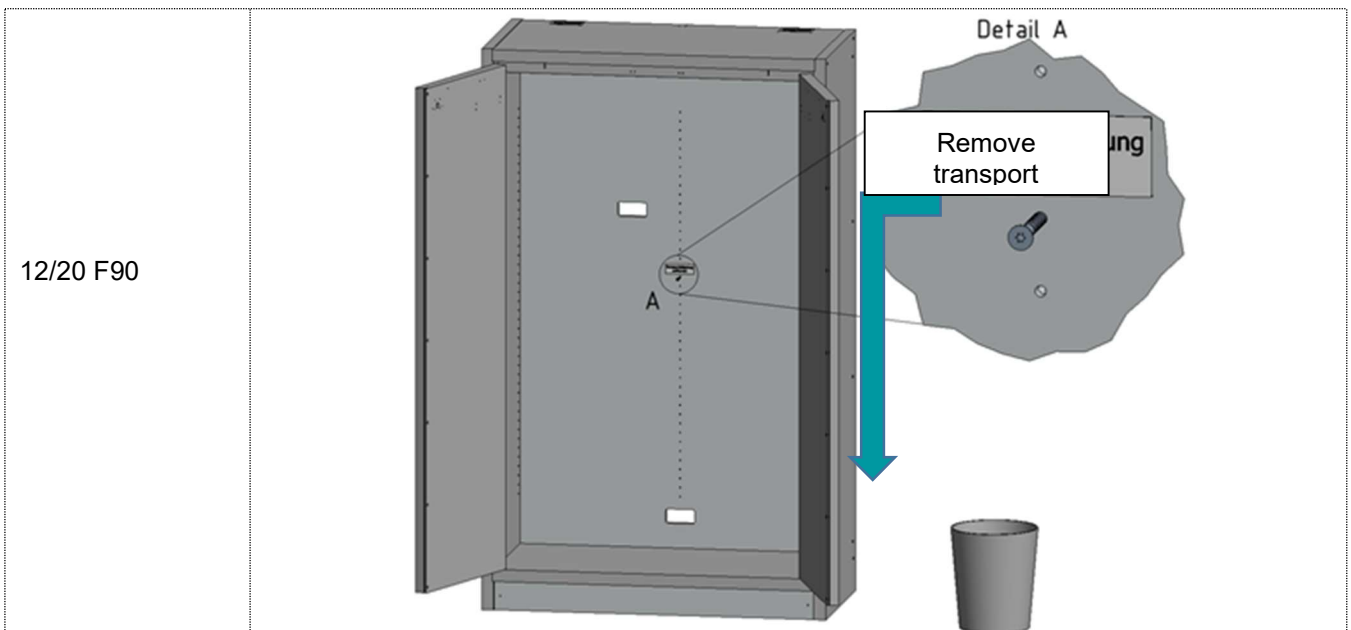Technical changes are reserved!

5 Closure types/Locks


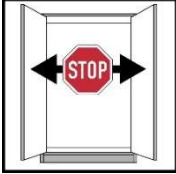
 Item	 Fig. 1	 Fig. 2	 Fig. 3
6/20 F30	✓	✓	✓
12/20 F30	✓	✓	✓
11/6 F30	✓	✓	✓
6/20 F90	✓	✓	✓
12/20 F90	✓	-	✓
11/6 F90	✓	-	✓

Fig. 1 The doors close automatically in case of fire.

Fig. 2 Doors with arrest system

Fig. 3 Lockable using a cylinder lock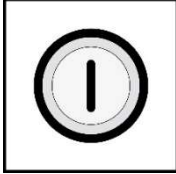



The products should in general always be kept closed when not in use. The products can be locked using a lock cylinder to protect against unauthorized access. It should be noted that the products have no emergency unlock, i.e. that persons can become locked in the cabinet unable to free themselves unaided!

6 Interior fittings

6.1 Base tray


Pursuant to EN 14470-1:

A base tray must be installed beneath the lowest setting level. The base tray must have a minimum collection volume of 10% of all of the containers stored in the cabinet or at least 110% of the volume of the largest individual container depending on which volume is greater.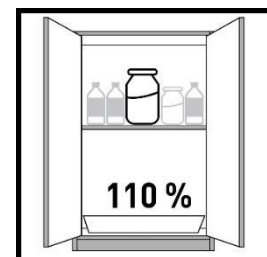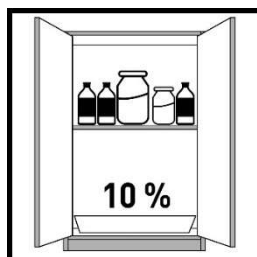

Attention:

Use of the base tray as a storage surface is permissible only in connection with a perforated plate insert.

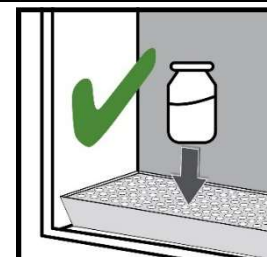Technical changes are reserved!

6.2 Tray bases

6.3 Pull-outs

Please obtain information from the enclosed operating manual here.

7 Storage


- Containers with aggressive chemicals (acids and bases) are to be accommodated in special acid/base cabinets or acid/base drawers as well as in safety cabinets with non-metallic interior fittings.
- Stored corrosive liquids can impair the function of the shut-off devices for supply and exhaust air.
- Each discontinued container reduces the minimum collection volume required under 6.1 relative to the total amount of stock.

Technical changes are reserved!


8 Earthing


An earthing connection is required depending on the product's customer-specific use. The decision concerning this is made via a risk analysis by the customer.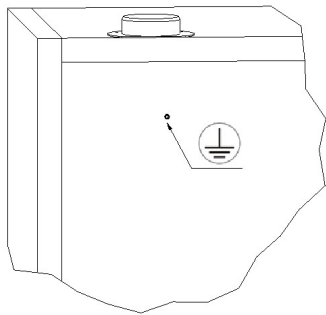


12/20 F90

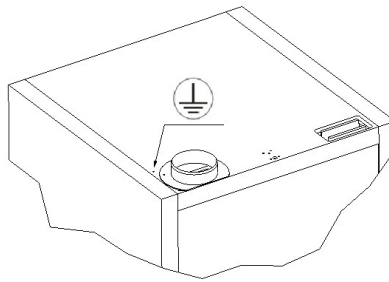
6/20 F90

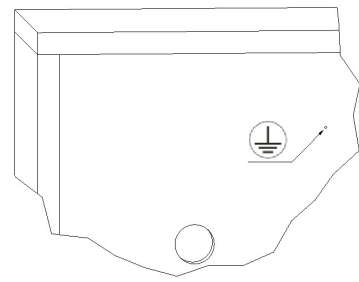
11/6 F90

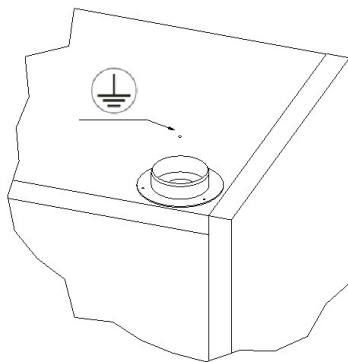12/20 F30

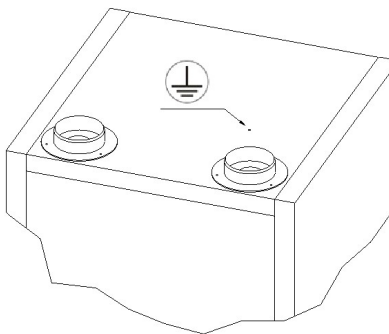6/20 F30

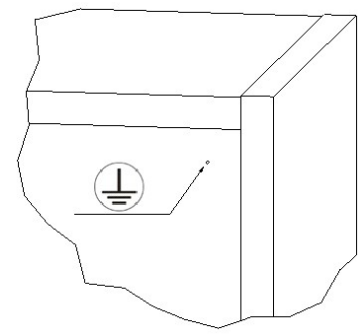
11/6 F30

Technical changes are reserved!

9 Ventilation

- All models can be connected via the air vent to a technical ventilation system. Obtain the air exchange prescribed for this, max. pressure loss, and exhaust guides from the respectively applicable national regulations.
- Care must basically be exercised to ensure that exhaust air is directed into a non-endangered area. PROTECTO recommends an external circulation cap if this cannot be guaranteed.

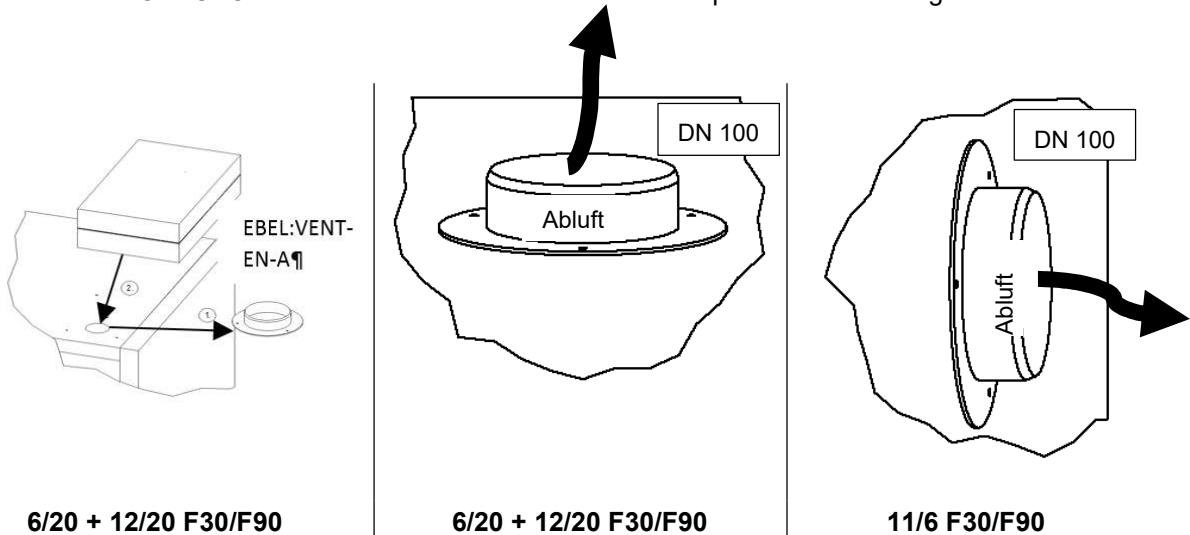Proper connection to a ventilation system must be inspected (e.g. using a flue tube). These products are generally also operated without technical air venting. However it must be noted here that the product's immediate vicinity might become an explosion-prone area. The user must attach a label to the product if it is to be used without an exhaust connection.

10 In case of fire

10.1 Opening the product after the fire

Combustible vapour-air mixtures may have formed depending on the fire's duration. Therefore remove all ignition sources in advance from a wide area around the product. Following a fire, the product may only be opened after six times the fire's duration. Only non-sparking tools may be used with extreme care for opening.


11 Disposal


The products can be disposed of correctly sorted. None of the components present needs to be sent to hazardous waste.

Technical changes are reserved!


12 Connection diagrams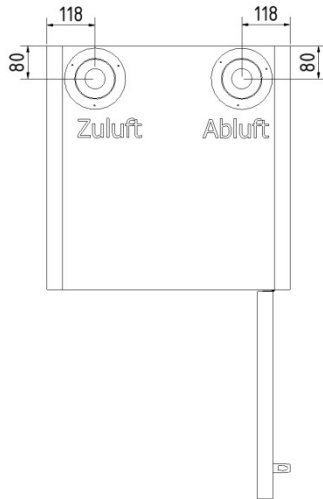


Type 6/20 F30 top view

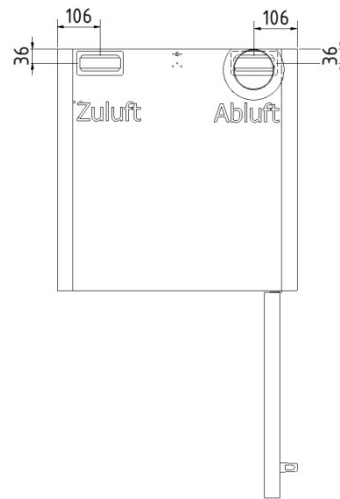
Type 6/20 F90 top view

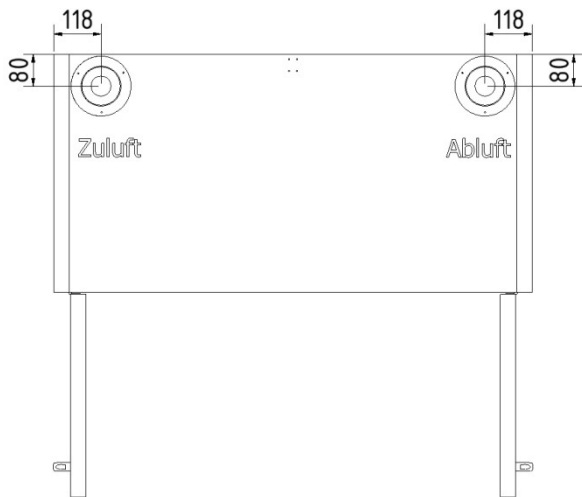
Type 12/20 F30 top view

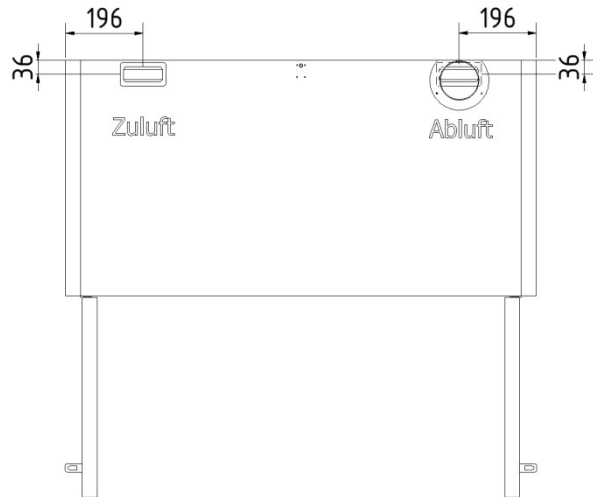Type 12/20 F90 top view

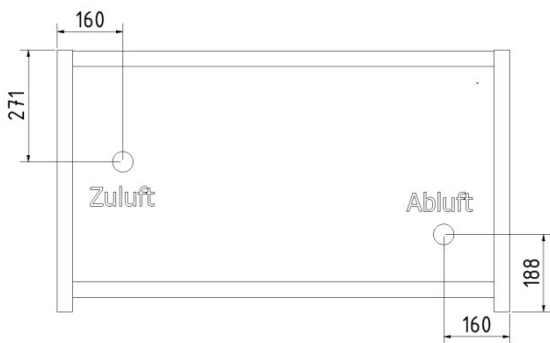Type 11/6 F30 rear view

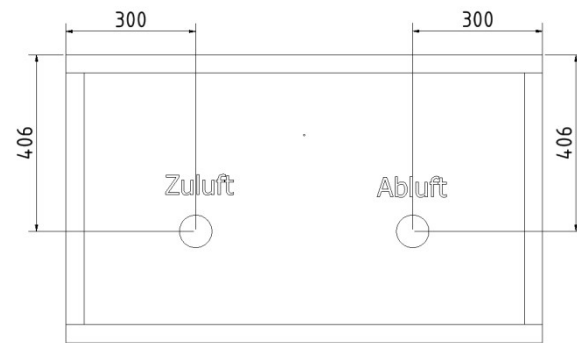Type 11/6 F90 rear view

Technical changes are reserved!

13 Optional accessories

13.1 General safety information

- Keep this Installation and Operating Manual adjacent to the cable / pipe bulkhead for each product.
- Affix any legally prescribed safety information at a location where it is conspicuous to users.
- Strictly observe:
 - All relevant VDE regulations
 - Fire authority regulations
 - General planning and building laws and regulations
 - Accident prevention and workplace regulations
 - Any instructions given by the technical supervisory service
- Pay attention to DIN 4102 regarding the assessment of fire protection and structural aspects.
- Work on electrical systems must be carried out with the power disconnected and performed by electricians.
- Never use a product that is not in perfect condition.
- Ensure that all required safety audits are carried out by an authorised expert.
- Damage or malfunctions due to incorrect transport or installation can only be reliably avoided if these activities are carried out by persons trained and authorised by us.
- Heed the technical information given in our catalogue and data sheet regarding the product concerned.
- Note that “hot spots” may form depending on the position of the fitted components. These components should therefore be distributed as evenly as possible. Check this if necessary by taking measurements inside the cabinet.
- Unauthorised persons must be denied access to the cable / pipe bulkhead.
- Incorrect installation could impair the protective function.
- Observe all information contained in this Installation and Operating Manual.
- Use only suitable, approved screws / dowels to fasten the cabinet to the building or structure.
- Make sure all screws / dowels are suitable for your particular building / structure.
- The doubler plates of the cable / pipe bulkhead must not be glued, painted or otherwise coated.
- There is not enough space in this manual to list every relevant regulation or directive. Please therefore make sure your installation is in conformity with all applicable regulations.

13.2 Intended use


- The product described in this Installation and Operating Manual was manufactured in accordance with the latest state of the art as well as generally recognised safety standards
- It may only be used:
 - For its intended purpose, and
 - If in perfect order and condition
- All other forms of use are considered to be contrary to the intended purpose. PROTECTO is not liable for any direct or consequential damages resulting from such use

13.3 Warranty, surrender of goods

- All installation instructions must be observed in order to guarantee optimal functioning of our products.
- PROTECTOPLUS cannot be held responsible for damages to other products which result from the installation or use of this product.
- The warranty which is given by PROTECTOPLUS is limited to products shipped by us.
- Design modifications carried out without prior consultation with PROTECTOPLUS result in voiding of the approval / warranty.
- Installation services must be guaranteed by the entity responsible for the installation.
- Warranty and liability claims for personal injury or property damage are excluded if they are the result of one or more of the following:
 - Improper installation, commissioning, operation or maintenance,
 - Failure to observe the information on transport, storage, operation and installation,
 - Inexpert repairs, or
 - Emergencies due to third party actions or force majeure.

13.4 Single-cable / single-pipe penetration

Scope of supply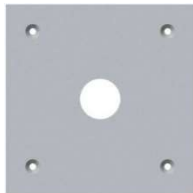

1x



4x

1x

Positioning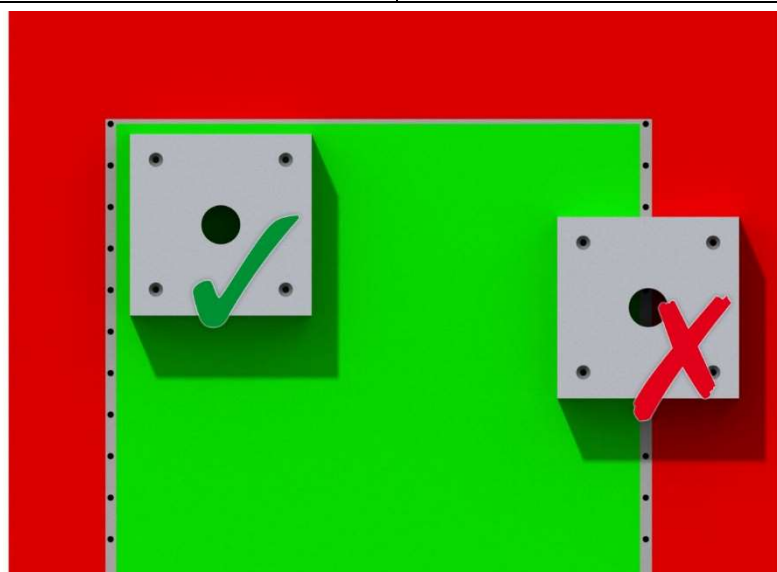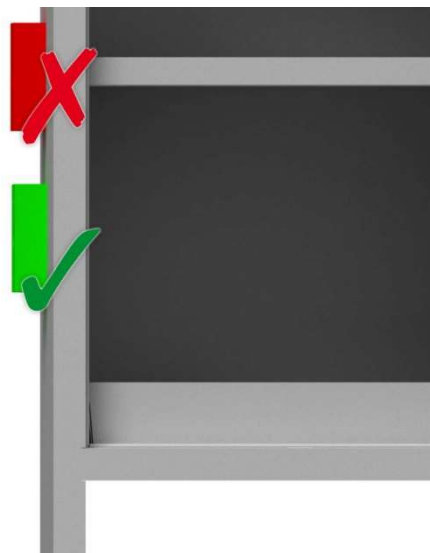

Installation procedure


1

2

3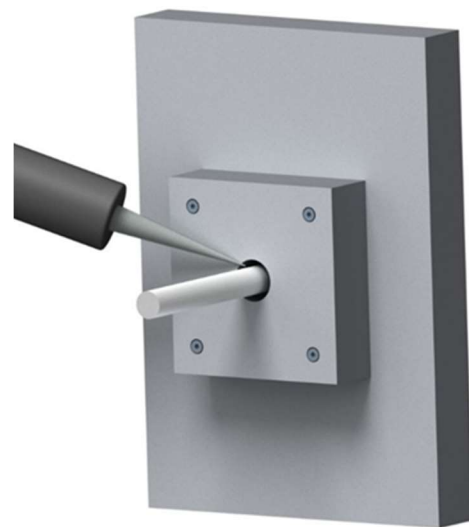

4

Technical changes are reserved!

13.5 Fire-resistant cable / pipe bulkhead

- The doubler plates are comprised of coated, non-metallic, plate-shaped components, which are painted light grey (similar to RAL 7035) on both the inside and outside. The edges are in a different colour.
- The fire insulation properties are achieved by processing fire protection panels to DIN 4102
- Provided they are installed in the proper manner, the cable / pipe bulkhead has a fire resistance rating of more than 30 or 90 minutes and is smoke-tight in the presence of (smouldering) cable fires.
- The information contained in this manual should be strictly observed to ensure that the product and the cable entry are handled correctly.
- Max. cable occupancy: 60%

Scope of supply

- 3x doubler plate 200 x 200 mm, with central through-hole Ø135 mm and countersunk fixing holes in the four corners
- 12x fixing / countersunk head screw 4.5 x 70 mm
- 2x firestop plug Ø 134 mm
- 1x firestop silicone cartridge
- 1x Installation and Operating Manual

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- The position of the cable / pipe bulkhead (on our products) should be
- determined on site, taking the dimensions of the doubler plate into account
- Drill a Ø 135 mm hole with a core drill (see Figure 1)
- The doubler plates can be used as a marking template

Figure 1

- Smooth the drilled hole with sandpaper (CAUTION: The edges are sharp)
- Mark and drill Ø 3.5 x 35 mm pilot holes on site for fixing the doubler plates (see Figure 2)

Figure 2

- Position the doubler plates exactly over the \varnothing 135 mm hole, then fasten them using the enclosed countersunk head screws after drilling the pilot holes in the wall
- The fixing holes in the doubler plate are offset, so that one plate can be screwed to another if the plates are turned 180°
- After fastening the doubler plates, insert the firestop plugs (one each on the inside and outside (see Figure 3)).

Figure 3

- Feed the cables through, then seal the openings with the silicone which is supplied with the product.

- F30 cable / pipe bulkheads: Two doubler plates plus the wall plate on which the bulkhead is installed -> wall thickness 120 mm
- F90 cable / pipe bulkheads: Three doubler plates plus the wall plate on which the bulkhead is installed -> wall thickness 150 mm
- Note that the required wall thickness also includes a layer of internal insulation (approx. 40 mm)

13.6 Recirculation fan, type SST-UL V4

13.6.1 Instructions for safe use

To ensure the safe operation of the system, the following basic requirements must be met:

- Temperature at place of installation: between +10°C and +40°C
- Maximum humidity at place of installation: 55%
- Limit value for dust content in ambient air: 10mg/m³

The SST-UL recirculation fan is used to extract and clean the exhaust air from hazardous material, safety, chemical and gas cylinder cabinets. The device ventilates the cabinet depending on the required air change rate, filters out hydrocarbons and indicates the current function by means of two lamps each (RED = Error, GREEN = Service). The concentration of hazardous substances in the exhaust air is shown in 25% steps. A concentration of 100% corresponds to the maximum load. The DSA (dynamic saturation adjustment) function corrects the switching point according to the ambient air quality. The device also has two potential-free outputs for these signals. It is suitable for extracting air from cabinets as defined in EN 14470-1 and DIN 12925 Part 2 and conforms to the requirements of TRGS 510 and GefStoffV. The user is responsible for ensuring safe operation of the system. The statutory limit values (TRGS 900 "Occupational exposure limit values" and TRGS 903 "Biological limit values") must be observed at the place of use.

PERSONAL INJURY AND MATERIAL DAMAGE

Improper use could result in broken impellers, fatigue failures or fires due to sparks, possibly leading to personal injury or material damage.

13.6.2 Specific hazards

The device conforms to 2014/34/EU (ATEX) (equipment group II, category 3 and temperature classification T4); it is exclusively intended for extracting air from Zone 2 or non-Ex zones and must be installed in a safe area.

Marking:   II 3/-G Ex ec IIC T4 Gc

All forms of use which deviate from or exceed the limits of use described above are considered to be contrary to the intended purpose. The manufacturer / supplier is not liable for any damages resulting from such use; the sole responsibility lies with the user.

No specific hazards for people or the environment arise provided this equipment is used as intended and only with original accessories, and provided this Installation and Operating Manual as well as the information on safe commissioning are observed.

13.6.3 Operating principle, safe installation, commissioning, adjustment, installation / removal, maintenance

The device is shipped ready to use and install with the exception of the control cables, which are not included in the scope of supply because they must be adapted to the specific conditions on site. You must check the electrical specifications (voltage, frequency) and compare them with the information on the rating plate prior to connecting the device to the mains.

The device may only be installed, operated and maintained by appropriately trained, instructed and authorised persons.

- Operating principle: The air containing hazardous substances is discharged from the hazardous material cabinet via a filter system comprised of a coarse filter, a fine filter and an activated carbon filter. 99.9% of all hydrocarbons are retained in this way by adsorption. This micro-controlled device monitors the volume flow and saturation of the activated carbon filter. The operational status is indicated by three LEDs (GREEN, YELLOW and RED) as well as by an audible signal. The exhaust function is interrupted as soon as the saturation LED shows "Error". This ensures that no hazardous substances are discharged from the cabinet into the workroom. The interior of the device is automatically flushed in the event of an error in order to reduce the concentration. If flushing does not cause the concentration to fall below the setpoint, the device goes into "alarm". Alarms can only be cancelled by authorised persons. Short concentration peaks (less than 9 seconds) are simply indicated (all saturation lamps on) without the device going into "error". The coarse and fine filters significantly extend the life of the activated carbon filter by retaining dust particles. The actual saturation up to the limit value is indicated by the three yellow LEDs in 25% steps. The quality of the ambient air is shown by another yellow LED. The brightness of this LED varies according to the amount of the hazardous substances.
- Installation location: the capacity of the recirculating air filter is intended for cleaning the air volume from the cabinet (cabinet volume x air turnover). Since the exchange air is taken from the ambient air via cabinet air intake, it must be taken into account that the concentration of pollutants is influenced by the composition of the ambient air. The total concentration of pollutants before the filter is therefore composed of the sum of the cabinet and ambient concentrations. When installing the cabinet, make sure that the ambient air is pollutant-neutral. The device provides an indication of the quality of the ambient air using the four-level


"DAS" brightness display (in the normal state, the LED does not light up). The brightness levels are visible in test mode - operation of the test button.

- Polluted air circulation can result in frequently occurring fault messages and can eventually lead to shutdowns due to increased concentration values. In order to ensure the normal operation of the system, the following installation situations should be avoided:
 - Proximity of cleaning facilities where solvent or cleaning agents are handled.
 - Proximity of storage of outgassing substances.
 - Proximity of workplaces that generate a lot of dust (sanding, polishing, sand blasting etc.)
 - Proximity of workplaces that generate a lot of moisture (washing facilities, cooking facilities, chemical/thermal surface treatment baths, dryers, etc.)
 - Proximity of nearby painting activity
 - Proximity of exhaust gas emissions (engine tests in repair shops, forklift traffic, incinerators etc.)
- Installation and commissioning: The device is attached to the cabinet by means of the connection fitting (accessory). The potential-free signal outputs are connected by the customer if necessary. The device goes into test mode as soon as it is connected to the mains supply with the mains plug. The LEDs, potential-free outputs and sensors are then checked one at a time.
- Function check: A function check should be carried out whenever the device is started up or as needed (see "Troubleshooting"). When the device is connected, or if you press the TEST button, all lamps light up for 3 seconds. Following this, the relays for the potential-free contacts are switched on and off for 3 seconds each. The flow and saturation sensors are checked in the final test sequence. The flow is checked by switching the technical ventilation system on and off (5 seconds on, then 5 seconds off). The LEDs for the flow change colour from RED to GREEN and the result of the test is indicated during the pauses (RED = Error, GREEN = OK). At the end of the test, the device automatically goes into normal mode. The exhaust air monitoring function can also be checked by briefly covering over the exhaust vent in order to throttle the volume flow.
- Adjustment: The exhaust air monitoring function is set to the required flow and saturation at the factory, so that no further adjustments are necessary when you start up the device for the first time.
- Repairs: Repair work may only be carried out by authorised persons.
- Maintenance: The device must be checked by an authorised person as and when necessary (once a year as a minimum). The filters must be replaced or cleaned at the same time. Replacement filters can be ordered from the supplier. We recommend stocking a set of replacement filters to avoid interrupting operation. The device records the operating time internally and outputs an audible signal (one 3-second beep every hour after 365 days) to indicate that maintenance is due.


13.6.4 Troubleshooting

Fault	Cause	Remedy
Device not working	No power Fuse defective Device defective	Check the mains voltage Replace the fuse of the mains connector Send in the device
Red flow lamp ON	Dust filter clogged Volume flow interrupted	Replace the coarse filter Replace the fine filter Remove the obstruction from the exhaust vent
Red saturation lamp ON, audible signal, air flow interrupted	Activated carbon filter saturated Concentration of hazardous substances too high Other hazardous substances contained in exhaust air	Replace the filter Check the contents of the cabinet and remove any leaked substances Start up the device again (authorised persons only) Contact the supplier


13.6.5 Specifications, characteristics

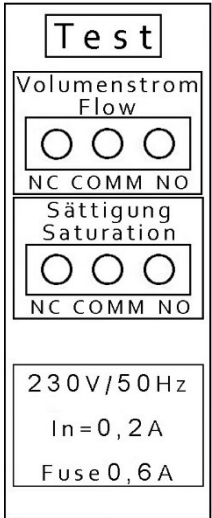Model	: SST-UL	
Ventilation system	: RCRv75/2m	
Flow	: 10 to 25 m ³ /h	
Total pressure	: Max. 250 Pa	
Sound level	: 38 dB(A) at a distance of 3 m	
Monitored parameters	: - Flow - Saturation of activated carbon filter	
LEDs	: ROT = Error, GREEN = Service 25 / 50 / 75% Ambient concentration	
Signal outputs	: - Potential-free for flow - Potential-free for saturation of activated carbon filter	
Contact power	: 230 V / 8 A	
Power supply	: 230 V / 50 Hz	
Current consumption	: Max. 0.5 A	
IP protection	: CE Ex II 3/-G Ex ec IIC T4 Gc PTB 21 ATEX D137	
Frequency	: 50 Hz	
Motor protection	: PTO 90°C	
Weight	: 20 kg	


Characteristic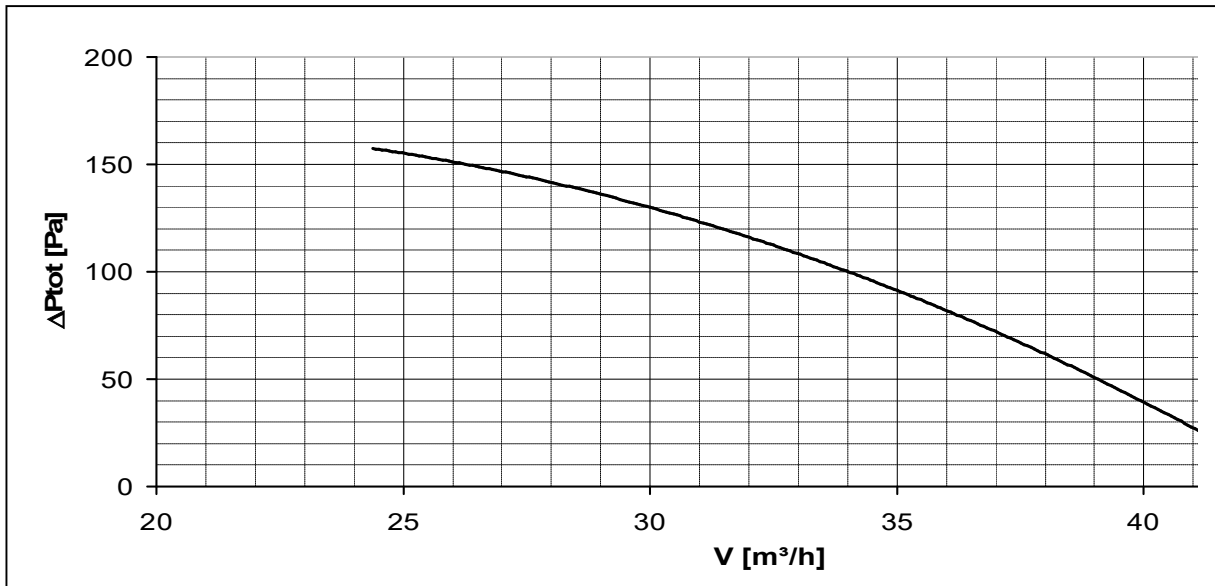

Dimensions

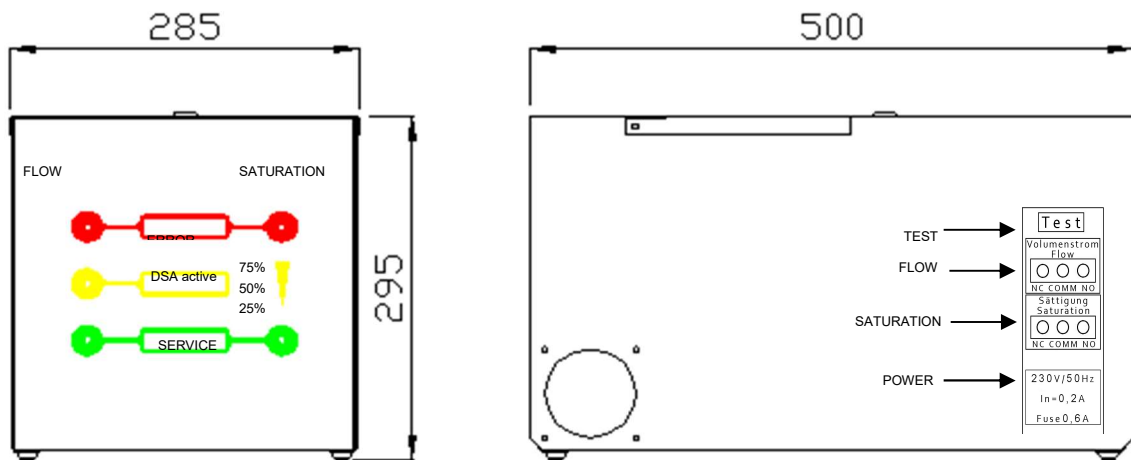13.6.6 Accessories shipped with the SST-UL

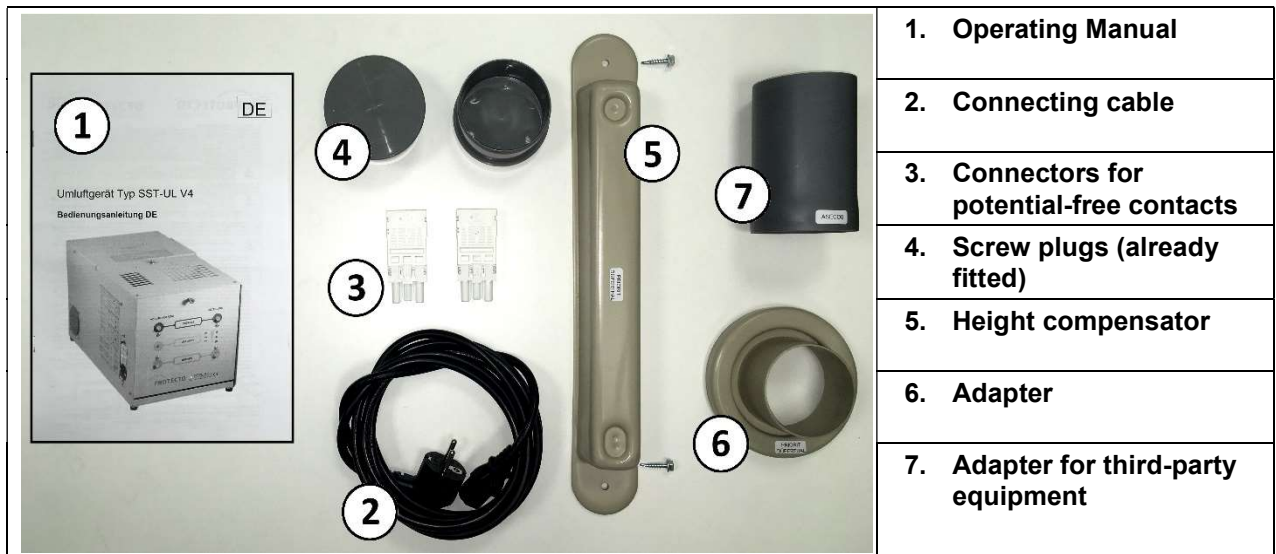13.6.7 Installing the SST-UL

- Insert the adapter
- Place the device on top of the cabinet with the adapter over the exhaust vent, then place the height adjuster underneath the device and align
- Fasten the height adjuster to the cabinet using the screws provided

13.7 Exhaust fan, type SST AL

13.7.1 Instructions for safe use

The **SST-AL** exhaust fan is used to extract hazardous substances from hazardous material, safety, chemical and gas cylinder cabinets as well as from acid and lye cabinets (aggressive and / or explosive media with a dust content of $< 5 \text{ g/m}^3$). Ambient temperatures from $+10^\circ\text{C}$ to $+40^\circ\text{C}$ are permitted. The device ventilates the cabinet depending on the required air change rate and indicates the current function by means of two lamps each (RED = Error, GREEN = Service). The device also has potential-free outputs for these signals. It is suitable for extracting air from cabinets as defined in EN 14470-1 and DIN 12925 Part 2.

PERSONAL INJURY AND MATERIAL DAMAGE

Improper use could result in broken impellers, fatigue failures or fires due to sparks, possibly leading to personal injury or material damage.

13.7.2 Specific hazards

The device conforms to 2014/34/EU (ATEX) (equipment group II, category 3 and temperature classification T4); it is exclusively intended for extracting air from Zone 2 or non-Ex zones and must be installed in a safe area. Marking: **CE Ex II 3/-G Ex ec IIC T4 Gc**

All forms of use which deviate from or exceed the limits of use described above are considered to be contrary to the intended purpose. The manufacturer / supplier is not liable for any damages resulting from such use; the sole responsibility lies with the user.

No specific hazards for people or the environment arise provided this equipment is used as intended and only with original accessories, and provided this Installation and Operating Manual as well as the information on safe commissioning are observed.

13.7.3 Safe installation, commissioning, adjustment, installation / removal, maintenance

The device is shipped ready to use and install with the exception of the control cable and exhaust duct, which are not included in the scope of supply because they must be adapted to the specific conditions on site. You must check the electrical specifications (voltage, frequency) as well as the resistance of the system and compare them with the information on the rating plate prior to connecting the device to the mains.

The device may only be installed, operated and maintained by appropriately trained, instructed and authorised persons.

- **Installation and commissioning:** The device is attached to the cabinet by means of the connection fitting (included). The signal output and the exhaust duct are connected. You must ensure that no foreign objects can get into the exhaust air monitoring system via the exhaust duct (protective grid). The device is ready to use as soon as it is connected to the mains supply with the mains plug. The green lamp indicates normal mode.
- **Function check:** A function check should be carried out prior to connecting the exhaust duct or as needed (see "Troubleshooting"). To do this, attach the device and connect it to the mains supply. The device is ready to use as soon as it is connected. The green lamp now lights up permanently. The device should go into error mode if you cover over the exhaust air outlet briefly in order to throttle the volume flow. Repeat this procedure several times. In order to test the monitoring function without removing the device, press the TEST button on the rear.
- **Adjustment:** The exhaust air monitoring function is set to the required flow at the factory, so that no further adjustments are necessary.
- **Installation / removal:** Repair work may only be carried out by authorised persons.
- **Maintenance:** The device requires no maintenance.


13.7.4 Troubleshooting

Fault	Cause	Remedy
Device not working	No power Device defective	Check the mains supply Send in the device
Red lamp permanently ON or blinking, function check OK	System resistance too high	Change the duct layout, reduce the air velocity in the duct (by increasing the cross-section)
Red lamp permanently ON, function check not OK	Ventilation system sticks, control loop defective	Send in the device

13.7.5 Specifications

Model	:	SST-AL
Ventilation system	:	RCRv75/2m
Ventil. housing material	:	Acid and lye-resistant plastic
Impeller material	:	Acid and lye-resistant plastic
Flow	:	50 to 200 m ³ /h
Total pressure	:	250 to 50 Pa
Speed	:	2400 to 2800 rpm
Sound level* Lp2A	:	35 dB(A)
LEDs	:	ROT = Error, GREEN = Service
Operating cycles	:	Min. 1,000,000
Signal output	:	Potential-free via DIN bush
Contact power	:	Max. 230 V / 8 A
Power supply	:	230 V +5/-10%
Current consumption	:	Max. 0.45 A
Frequency	:	50 Hz
Motor protection	:	PTO 90°C
IP protection	:	CE Ex II 3/-G Ex ec IIC T4 Gc PTB 21 ATEX D137
Weight	:	8 kg

Signal output assignment
Max. 230 V / 8 A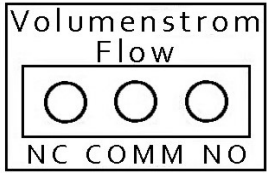


Volumenstrom
Flow


NC COMM NO

* At a distance of 3 m in accordance with DIN 45635, with air connections on the suction and discharge sides

Characteristic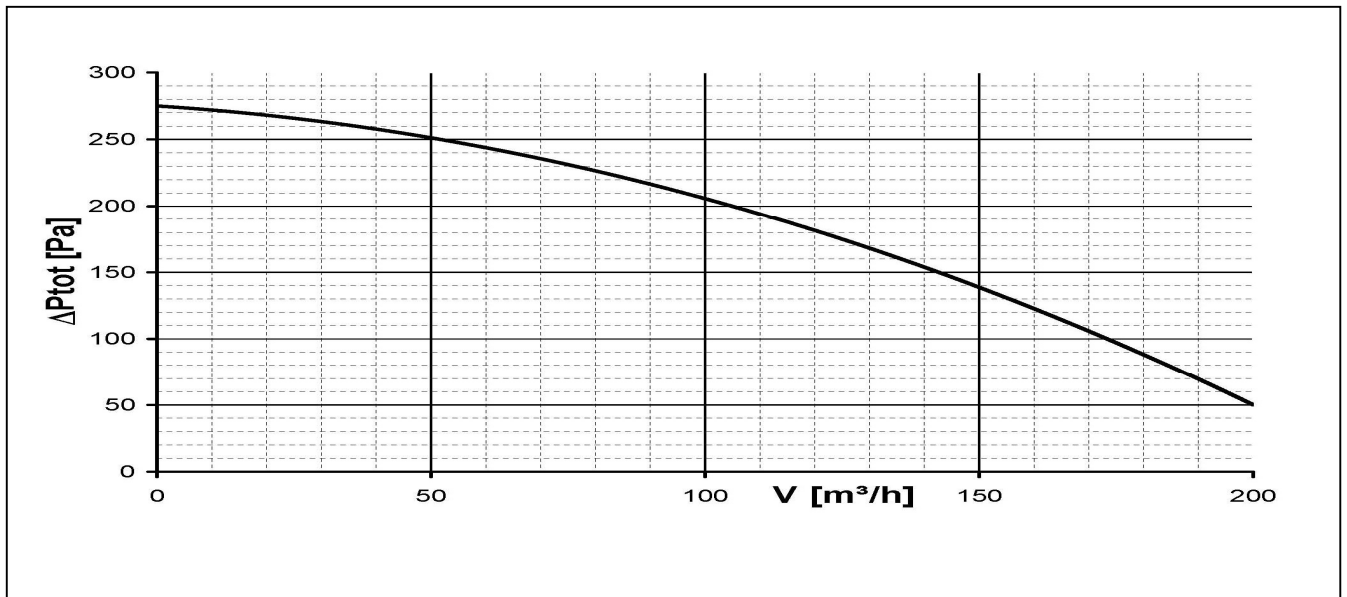

Dimensions

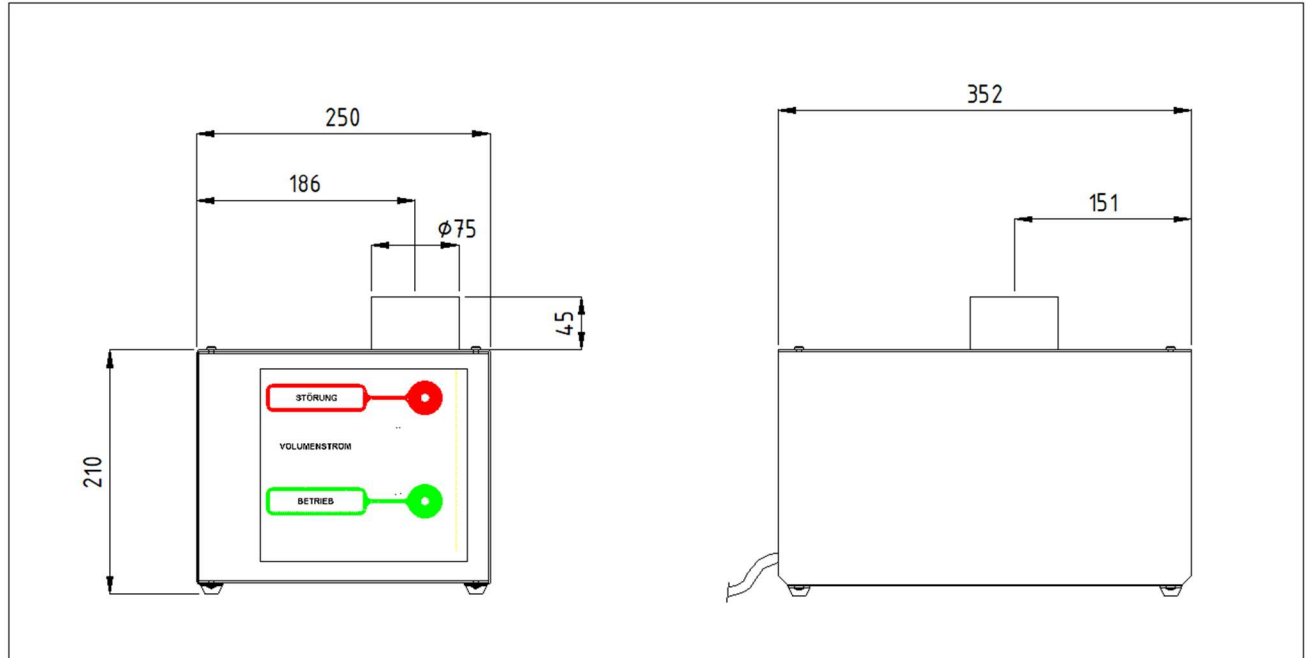13.7.6 Accessories shipped with the SST-AL

	1. Operating Manual
	2. Connectors for potential-free contacts
	3. Connection fitting for all cabinet types, already fitted
	4. Connection fitting (type SST-D only)
	5. Adapter for third-party equipment
	6. Adapter for SST-P

13.7.7 Installing the SST-AL for cabinets

- Insert the adapter (6) at the free end of the connection fitting and fasten it with the hose clamp provided
- After preparing the adapter in this way, insert it into the device and the exhaust air outlet of the cabinet and fasten it to the cabinet outlet using the hose clamp provided. Align the device.

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MPA Dresden GmbH